
MASTERflex®
(US & Canada only) Toll Free 1-800-MASTERFLEX • 1-800-637-3739
(Outside US & Canada) 1-847-381-7050
www.masterflex.com • techinfo@masterflex.com
*EN809 Manufactured by:
Cole-Parmer Instrument Company
28W092 Commercial Avenue
Barrington, IL 60010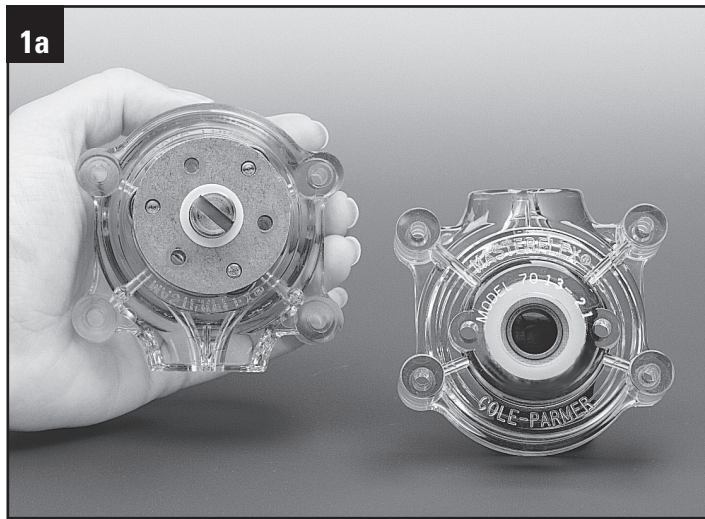

1a
ENGLISH

1. Single Pump Head Loading

⚠ WARNING: PRODUCT USE LIMITATION

These products are not designed for, nor intended for use in patient connected applications; including, but not limited to, medical and dental use, and accordingly have not been submitted for FDA approval.


Note: Use only MASTERFLEX precision tubing with MASTERFLEX pumps to ensure optimum performance. Use of other tubing may void applicable warranties.

Contents: One pump head, one 15 in (38-cm) length of silicone tubing, one mounting hardware package, manual and tubing loading key.

Supplied tubing loading key required for assembly.

- Separate the end bells (the pump head halves). Hold the end bell containing the rotor as shown with the tubing retainer grooves pointing down.
- Place tubing in the right groove and against the first two rollers. Hold tubing with thumb. Near groove, insert smaller prong of loading key between the top of the rotor and tubing. Push key in as far as possible.
- Push down and turn key counterclockwise (ccw) completely around the rotor. The key will push the tubing uniformly into the end bell assembly. Hold the second end of tubing. Remove the key.
- Position the other end bell on top and press the end bells together. Be careful not to pinch the tubing. If end bells do not snap tightly together, reload tubing. If necessary, turn key in slot on rotor shaft to adjust tubing (as in Step e).
- With key in slot on rotor shaft, turn key to align tang on rotor shaft with slot in motor drive shaft. Point tubing retainer grooves up. Shift the pump head slightly until it snaps on the alignment pins (if present). Secure with four provided screws. Tighten with fingers only.

PUMP FOR LIQUIDS ORIGINAL INSTRUCTIONS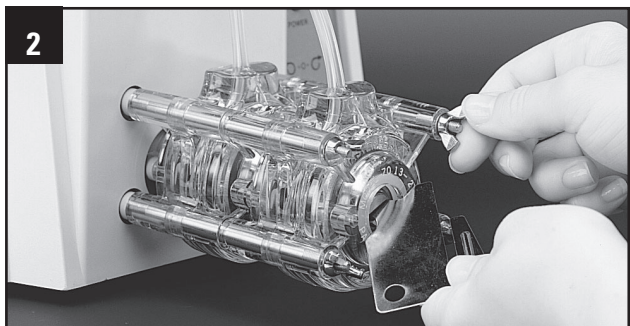

2. Multi-channel Mounting


Flat bladed Screwdriver required for mounting.

Tubing loading key required for mounting.

Note: Order special mounting hardware for multi-channel pumping, see "3. Replacement Parts and Accessories".

- Load the pump heads with tubing.
- Install the four correct length mounting screws into the drive.
- Slide the first pump head onto the mounting screws.
- Place key in slot on rotor shaft. Twist to align tang on rotor shaft with slot in motor drive shaft. Shift the pump housing around until it drops over the alignment pins (if present).
- Repeat for each additional pump head, aligning pump head tang with slot on previously mounted pump head.
- Slide the four flat washers onto screws and secure with the four wing nuts. Tighten with fingers only.
- A support bracket is supplied with 3- and 4-channel mounting hardware for additional support. Mount over bottom two screws. Insert one of three different adjustment screws depending upon drive height.

Trademarks bearing the © symbol in this publication are registered in the U.S. and in other countries.

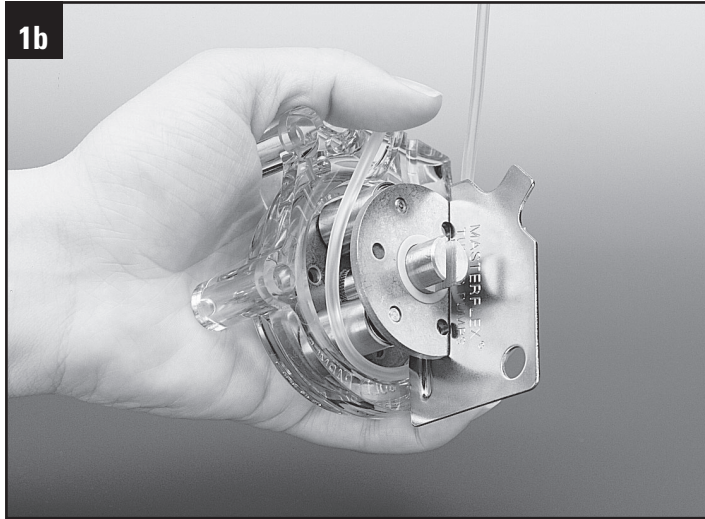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
A-1299-0108B Edition 23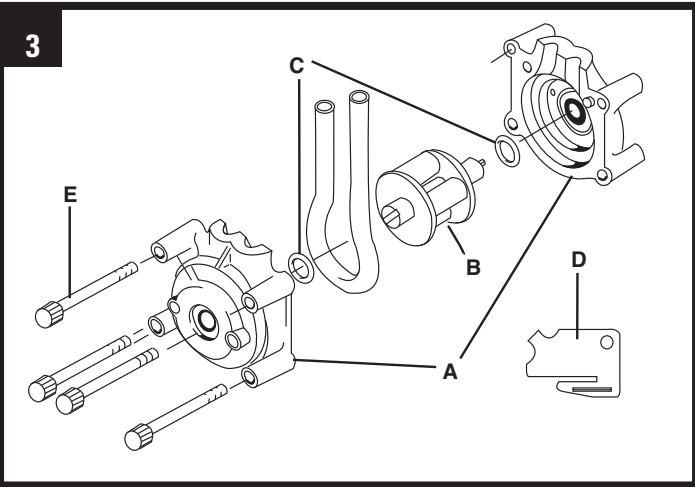

ENGLISH

3. Replacement Parts and Accessories

A. End Bells (order two end bells for a complete head assembly).

Pump Head number	PC Order number	Pump Head number	PPS Order number
07013-00, -20	07013-81	—	—
07013-10, -21	07013-91	—	07013-92
07014-00, -20	07014-81	—	—
07014-10, -21	07014-91	—	07014-92
07015-00, -20	07015-81	—	—
07015-10, -21	07015-91	—	07015-92
07016-00, -20	07016-81	—	—
07016-10, -21	07016-91	—	07016-92
07017-00, -20	07017-81	—	—
07017-10, -21	07017-91	—	07017-92
07018-00, -20	07018-81	—	—
07018-10, -21	07018-91	—	07018-92
07024-00, -20	07024-81	—	—
07024-10, -21	07024-91	—	07024-92
07035-02, -20	07035-81	—	—
07035-12, -21	07035-91	—	—

Pump Head number	Pump Head suffix	Order number
07013, 07014, 07016, 07017, 07018	-00 -10, -50 -20 -21, -52	07013-75 07013-76 07013-80 07013-95
07015, 07024, 07035	-00, -02 -10, -50, -12 -20 -21, -52	07015-75 07015-76 07015-80 07015-90

C. 07021-04 Thrust washers. Pack of 10.

D. 07013-90 Tubing loading key.

E. Mounting hardware for standard pump heads.

Set contains four #8-32 screws, four washers, and four wing nuts.

Number of heads to be mounted	Stainless steel order number
1	07013-04
2	07013-05
3	07013-08
4	07013-09

4. Specifications		
	Thin wall*	Thick wall*
Maximum continuous discharge pressure—psi (bar):	20 (1.4)	25 (1.7)
Maximum intermittent discharge pressure—psi (bar):	35 (2.4)	40 (2.7)
Maximum vacuum:	660 (510 ¹) mm Hg; 26 (20 ¹) in Hg	
Maximum suction lift:	8.8 (6.7 ¹) m H ₂ O; 29 (22 ¹) ft H ₂ O	
Number of rollers:	3	
Occlusion:	Standard fixed	
Maximum pump speed (rpm):	600	
Nominal torque load:	6.5 kg-cm (90 oz-in)	
Housing materials:	Polycarbonate (PC) all models, or Polyphenylene sulfide (PPS) all models except 07035	
Roller/rotor materials:	Cold-rolled steel (CRS) or Stainless steel (SS)	
Operating temperature [†] :	0 to 40°C (32 to 104°F)	
Noise Level:	< 70 dBa @ 1 meter	
Compliance (for CE mark):	EN809 (EU Machine Directive)	

* Thin wall: tubing L/S 13, 14, 16, 17, 18. Thick-wall: tubing 15, 24, 35.

[†]With tubing 17 & 18.

‡ Use in this temperature range for continuous duty operation with no decrease in performance or product life. Pump heads will work outside this range with some possible reductions in performance or product life.

5. Warranty and Return of Items

Warranty

Use only MASTERFLEX precision tubing with MASTERFLEX pumps to ensure optimum performance. Use of other tubing may void applicable warranties.

The Manufacturer warrants this product to be free from significant deviations from published specifications. If repair or adjustment is necessary within the warranty period, the problem will be corrected at no charge if it is not due to misuse or abuse on your part, as determined by the Manufacturer. Repair costs outside the warranty period, or those resulting from product misuse or abuse, may be invoiced to you.

The warranty period for this product is two (2) years from date of purchase.

Product Return

To limit charges and delays, contact the seller or Manufacturer for authorization and shipping instructions before returning the product, either within or outside of the warranty period. When returning the product, please state the reason for the return. For your protection, pack the product carefully and insure it against possible damage or loss. Any damages resulting from improper packaging are your responsibility.

Technical Assistance

If you have any questions about the use of this product, contact the Manufacturer or authorized seller.

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